

#### Terrific Validation, Tech Support and Prompt Delivery

#### Data Sheet

#### Protease Inhibitor Cocktail (EDTA-Free, 100X in DMSO)

A blend of pan-protease inhibitors for protection of protein integrity.

Catalog No.	Size	Price
K4000	1 mL	USD 37
	100 μL × 30	USD 147
	1 mL × 10	USD 270
	1 mL × 100	USD 1700

## **Product Overview**

Protease Inhibitor Cocktail is a blend of pan-protease inhibitors for protection of protein integrity. Endogenous proteins are produced and removed in a balanced state, so their cellular levels are generally stable under stable environmental conditions. However, protein production is greatly halted and degradation is enhanced when cells are studied in vitro. To prevent the degradation of proteins under such conditions, one can utilize a cocktail of small molecule inhibitors to block the action of proteases.

## **Price Comparison**

Brand	100mL(1x)	1000mL (1x)
Selleck	37	270
Sigma	63	450
PIERCE	68	483
Roche	84	456

### **Contents**

Ingredient	100x Conc.	Target
AEBSF	104 mM	Serine proteases
Aprotinin	80 μΜ	Serine proteases
Bestatin	4 mM	Aminopeptidases
E-64	1.4 mM	Cysteine proteases
Leupeptin	2 mM	Serine and cysteine proteases
Pepstatin A	1.5mM	Acid proteases

You can order by phone, email or fax

Tel: +1-832-582-8158 Fax: +1-713-796-9816

Email: sales@selleckchem.com

Last updated on December 4, 2013

1/3

### **Product Details**

**Intro:** A blend of pan-protease inhibitors for protection of protein integrity.

Usage: Thaw at room temperature, add at 1:100 (v/v) dilution in solutions (such as cell lysate and

tissue extract), before Western Blot analysis, Co-IP, pull-down, IF, IHC, kinase assay and

etc...


Storage: -20°C 12 months.

Shipping: Blue ice

Pre-caution: Some DUB proteases (one example is ATAXIN-3) cannot be suppressed by traditional

protease inhibitors, such as E-64, AEBSF, bestatin, leupeptin and aprotinin. (Reference: Neil D. Rawlings, Guy Salvesen et al. Handbook of Proteolytic Enzymes, Vol.1, 2012).

### **Customer Reviews**

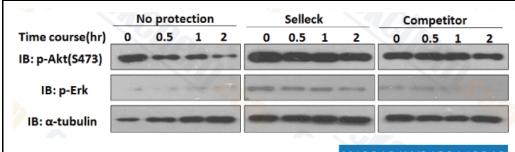

K400101W0120140313

Cell Lines 293T


Condition 0, 0.5, 1, 2hrs at room temperature

293T cells were harvested with three different lysis buffers, left at room temperature for indicated time, and then SDS-loading buffer was added into the lysate for Western Blot.

Results No Protection : EBC no phosphatase inhibitors

Selleck: EBC with protease & phosphatase inhibitor from Selleck (K4000,K5000)

Competitor : EBC with protease & phosphatase inhibitor from one competitor

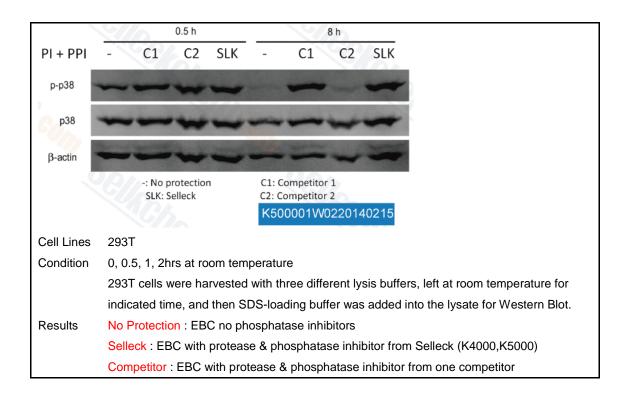

# **Related Cocktails**

### Phosphatase Inhibitor Cocktail (Full spectrum) K5000

## Quality guaranteed and expert technical support

Your money back if it does not work as we say on the datasheet.

For full details of the promise, please visit us online:

http://www.selleckchem.com/return-policy.html

You can order by phone, email or fax

Tel: +1-832-582-8158 Fax: +1-713-796-9816

Email:sales@selleckchem.com

Last updated on December 4, 2013